# Anti-Human P-cadherin, Monoclonal (Clone NCC-CAD-299)

**Code No. M127**

**Size : 0.1 mg Mouse Ig**

**Subclass : IgG1**

- \* 2 years from date of receipt under proper storage conditions.

#### **Source :**

Monoclonal antibody was obtained by fusing the mouse myeloma cell-line P3-X63-Ag8-U1 with spleen cells of BALB/c mouse after immunization with human epidermal carcinoma cell A431.<sup>1)</sup> The monoclonal antibody was harvested from ascitic fluid.

#### **Purification :**

Antibody was purified by column chromatography, dissolved in 10 mM PBS, pH 7.4, containing 1.0% bovine serum albumin, and then lyophilized.

The lyophilized antibody does not contain preservative.

**Form :** Lyophilized

#### **Reconstitution :**

Dissolve the lyophilized antibody in 50  $\mu$ l of distilled water (final concentration: 2.0 mg/ml). This solution can be used as a stock solution. If dilution is necessary for your application, dilute the stock solution with the following Dilution solution just prior to use. When the entire amount of antibody is to be used over a short time period, it may be dissolved directly in 500  $\mu$ l or more of the Dilution solution.

Note (1) : Be sure to store the antibody at a minimum concentration of 2.0 mg/ml. A lower antibody concentration may result in decreased stability.

Note (2) : Reconstituted antibody solution should contain 0.1% sodium azide as a preservative when stored at 4°C.

**Dilution solution :** 20 mM TBS (pH 7.5) \* 1  
10 mM CaCl<sub>2</sub>  
1.0% BSA  
(0.1% NaN<sub>3</sub>) \* 2

\* 1 TBS should be used for dilution, as PBS causes precipitation.

\* 2 When stored at 4°C, 0.1% sodium azide should be added as a preservative.

#### **Specificity :**

- This antibody specifically reacts with human P-cadherin.<sup>2),3)</sup>
- This antibody inhibits P-cadherin dependent cell-cell contact.<sup>2),3)</sup>

#### **Working concentration :**

- 100  $\mu$ g/ml  
(For inhibition assay of P-cadherin dependent cell-cell contact)
- 5 - 10  $\mu$ g/ml  
(For Western blotting under non-reducing condition by color detection)
- 5 - 10  $\mu$ g/ml  
(For immunohistochemical staining on frozen sections, by color detection)

#### **Storage :** 4°C

This product does not contain preservative.

The stock solution (2.0 mg/ml) should be stored in aliquots at -20°C for 1 year, or should be stored at 4°C for 6 months after adding 0.1% sodium azide. Avoid repeated freeze-thaw cycles. Diluted antibody should not be stored.

#### **Application :**

This product is sodium azide free and can be directly used for cell experiment.

- Immunohistochemical detection of P-cadherin on frozen tissue sections<sup>2)-6)</sup>
- Western blot analysis under non-reducing and non-heating condition (can not used for reducing condition)
- Analysis of P-cadherin dependent cell-cell contact mechanism

#### **References :**

- 1) Fabrictant R N, DeLarco J E, and Todaro G J. *Proc Natl Acad Sci USA.* (1977) **74:** 565-569.
- 2) Shimoyama Y, Hirohashi S, Hirano S, Noguchi M, Shimosato Y, Takeichi M, and Abe O. *Cancer Res.* (1989) **49:** 2128-2133.
- 3) Shimoyama Y, Yoshida T, Terada M, Shimosato Y, Abe O, and Hirohashi S. *J Cell Biol.* (1989) **109:** 1787-1794.
- 4) Shimoyama Y and Hirohashi S. *Cancer Res.* (1991) **51:** 2185-2192.
- 5) Yasui W, Sano T, Nishimura K, Kitadai Y, Ji Z-Q, Yokozaki H, Ito H, and Tahara E. *Int J Cancer.* (1993) **54:** 49-52.
- 6) Rasbridge S A, Gillett C E, Sampson S A, Walsh F S, and Millis R R. *J Pathol.* (1993) **169:** 245-250.

#### **Note :**

Ca<sup>2+</sup> should be contained in the buffers to stabilize cadherin antigen.

#### **Note**

This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc. Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from Takara Bio Inc.

If you require licenses for other use, please contact us by phone at +81 77 565 6973 or from our website at [www.takara-bio.com](http://www.takara-bio.com). Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements.

All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.

# Anti-Human P-cadherin, Monoclonal (Clone NCC-CAD-299)

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